Regulating transformer
12V/100VA

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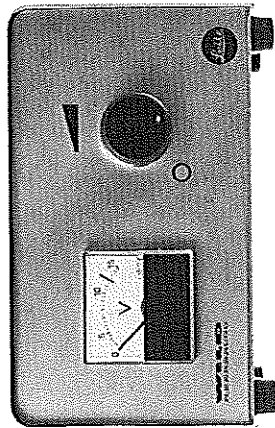
P6.2

Regulierttransformer
12V/100VA

Transformador
regulable 12V/100VA


10.156

WILD MTR27

Instructions for use
Mode d'emploi
Gebrauchsanweisung
Instrucciones para el uso

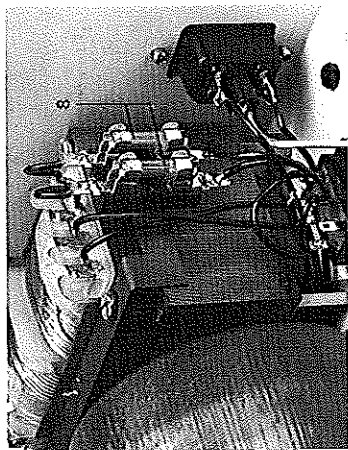WILD
HEERBRUGG

Echange des fusibles

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Changing the fuse
Disconnect the mains (line) cable, remove the two screws (7) from the underside of the transformer, and lift off the transformer housing. Replace the fuses (8) in the fuse holders. For all voltages from 110-240V, use 800 mA fuses. Replace and secure the housing.
Fuses: 2 X 800 mA, slow, 5 X 20 mm.

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

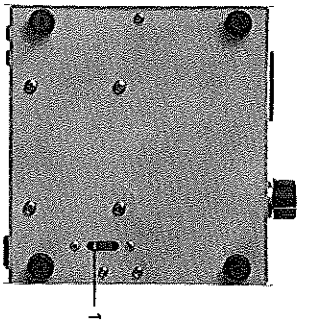
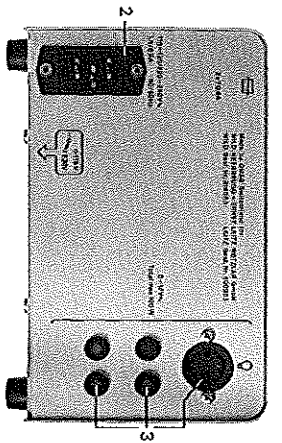
Setting the mains (line) voltage
Set the voltage selector (1) on the underside of the transformer to either 115V (for 110–120V) or 220V (for 220–240V). Plug the mains (line) cable into the socket (2) and connect it to the electricity supply.

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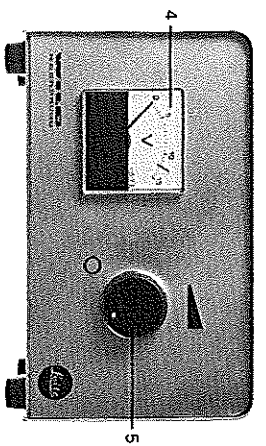Setting the click stop to the limiting voltage
Turn the regulating knob (5) clockwise to switch on, and then further until it engages the click stop. Slacken the screw (6) on the metal ring beneath the regulating knob, and turn the regulating knob (5) fully back (anti-clockwise).
Connect the 12V lamp to one of the three sockets (3). The total loading of these three sockets must not exceed 100W. Turn the regulating knob (5) until the voltmeter (4) reads 12V (red line). Retighten the screw (6) on the metal ring beneath the regulating transformer. Because the click stop is now at 12V, an accidental overloading of the lamp is avoided.
By the same procedure, the click stop can be set to 6V in order to run a 6V lamp.

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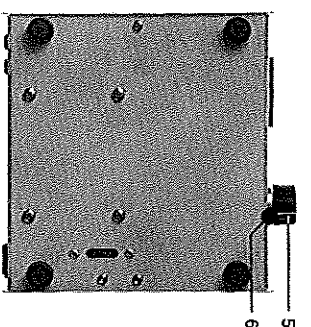New Transformer 12 V/100 VA (319 553)

Mains (line) voltage: Push voltage selector (S) for local mains (line) voltage.

Fuse: Through a corresponding circuit the transformer contains 2 identical fuses of 800 mA each for all mains (line) voltages from 110-250 V. Access to the fuses is gained by loosening the two cross-headed screws (indicated on the photograph by arrows).

Load: Since the transformer is laid out for 100 VA, the following lamps may be used with it:

Caution! Either 1 x 12 V/100 W (e.g., transmitted light, bright field-dark field) or 2 x 6 V/15 W low voltage lamps. Inspection of voltmeter: Do not greatly exceed 6 V voltage! Or 2 x 12 V/20 W (e.g., Epimakroscope M450).

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