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## Declaration of conformity C€

We the manufacturer Velp Scientifica s.r.l.  
Address Via Stazione, 16  
20865 USMATE (MB)  
Italy

under our responsibility declare that the product is manufactured in conformity with the following standards:

EN 61010-1 (2001)  
EN 61326-1 (2006)  
2002/95/EC (RoHS)  
2002/96/EC (WEEE)

and satisfies the essential requirements of the following directives:

Machinery directive 2006/42/EC  
Low voltage directive 2006/95/EC  
Electromagnetic compatibility directive 2004/108/EC  
plus modifications and that the documents listed in annex I are available at Velp's offices as foreseen by the machinery directive.

Velp Scientifica srl  
Via Stazione, 16  
20865 Usmate (MB) Italy  
Tel. +39 039 62 88 11  
Fax +39 039 62 88 120

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*Instruction for use: Laboratory stirrers type BS*

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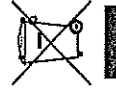
### IMPORTANT:

READ THE INFORMATION CONTAINED IN THE PRESENT MANUAL BEFORE USING THE UNIT.  
THE MANUFACTURER DOES NOT ACKNOWLEDGE ANY RESPONSIBILITY FOR AN IMPROPER USE OF THE EQUIPMENT, NOT RESPONDING TO DIRECTIONS FOR USE.

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
P 101.37

<b>Technical features:</b>		<b>BS</b>
Electric supply code for	230V/50-60 Hz	F20100151
Electric supply code for	115V/50-60 Hz	F20110151
Dimensions (BxHxP)	(mm)	80x215x196
Weight	(kg)	2,8
Power supply	(W)	60
Stirring speed	(rpm)	50 + 2.000
Speed display	(rpm)	NO
Resolution	(rpm)	-
Allowed room temperature	(°C)	0 + 40
Speed type	(rpm)	PWM
Safety against overheating		YES
Safety against overloading		YES
Safety display		-
Maximum stirrable volume (H <sub>2</sub> O)	(l)	25
Chuck with through hole Ø	(mm)	1 + 10
Admitted operation		Continuous

**Electric scheme:**

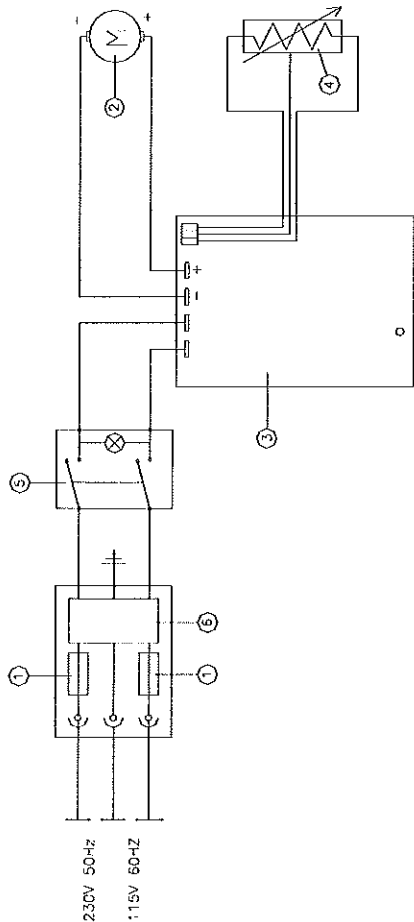- 1) Retarded fuse 5x20 (3, 15 A)
- 2) Motor
- 3) Electric board
- 4) Speed regulating potentiometer
- 5) General lighted switch
- 6) Filter

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